

Application Memo

Industry	Food
Instrument	Automatic potentiometric titrator
Measurement method	Acid-base titration
Standards	GB 1886.10 mmmø 0

In this memo, the concentration of the prepared 2w/w% acetic acid sample is tested. The diluted sample is titrated with 0.1mol/L sodium hydroxide up to pH8.3. The acetic acid concentration is calculated from the titration volume of sodium hydroxide.

Main unit	Automatic potentiometric titrator (preamplifier STD)
Electrode	Combined glass electrode Temmmmmmmø lufne